



System Requirements - Valab® Server V15.01



Characteristic	Possible values
Usage	Operation 24/7 (dedicated machine recommended, hibernation and network card sleep mode disabled). It communicates with your LIS through a TCP / IP network connection. Your LIS must have an option allowing Valab® connection and integration.
OS ⁽¹⁾	<ul style="list-style-type: none">• Windows Server 2019® x64 Standard Edition with Desktop Experience (graphical interface)• Windows Server 2016® x64 Standard Edition with Desktop Experience (graphical interface)• Windows Server 2012® R2 x64• Windows 10® x86 and x64 Pro edition
Processor	4-core - 64 bits - 2 Ghz
RAM	4 Gbytes
Disk type	Mechanical hard disk > 10 000 rpm or SSD
Typical recommended disk space	Space needed for the OS + : <ul style="list-style-type: none">- licence 1200 reports / day = 50 Gbytes- licence 2400 reports / day = 75 Gbytes- licence 4000 reports / day = 110 Gbytes- licence 6500 reports / day = 170 Gbytes- licence 10 000 reports / day = 250 Gbytes- licence 15 000 reports / day = 440 Gbytes
.NET Framework	Minimum 4.5.1 version
DB	No database engine installed
FireWall - Network	<ul style="list-style-type: none">• Fixed IP mandatory• Authorise incoming TCP protocol connections on the port(s) used to communicate with the LIS (by default port 7200 if only one LIS connection)• Authorise incoming TCP and UDP protocol connections on the ports used to communicate with the Valab® services of the Valab® Server and with the Valab® clients (by default 8080, 8089, 8090 and 2110)
Access	<ul style="list-style-type: none">• Local administrator account dedicated to the VALAB company• Telemaintenance access (VPN + RDP, remote control software,...)• Internet access to the "valab.com" domain
Backup / Security	<ul style="list-style-type: none">• Policy and implementation to be determined by the IS Manager of the medical laboratory• We recommend that you encrypt the connection between your LIS and Valab, in particular in the case of remote hosting of these systems.



System Requirements - Valab® Client V15.01

Characteristic	Possible values
Usage	Thick client for remote access to the Valab® server
OS ⁽¹⁾	<ul style="list-style-type: none">• Windows 10®• Windows Server 2019® x64• Windows Server 2016® x64• Windows Server 2012® R2 x64
Processor	dual core - 64 bits - 2 Ghz
RAM	4 Gbytes
Disk space	Space needed for the OS + 20 Gbytes
Graphics card	1366 * 768 minimum recommended
.NET Framework	Minimum 4.5.1 version
FireWall - Network	Open ports 2110, 8089, 8090 and 8080 to the Valab® server
Access	Read / write access to the installation directory of the Valab® client for the user

⁽¹⁾ Subject to effective application of the latest security updates provided by Microsoft® for the OS
On condition that the OS is still maintained by Microsoft®