

Valab® 16.02

Release Notes

The VALAB company is certified ISO 9001 by:



- Contents -

1	Object	3
2	Evolutions	4
2.1	LDAP connection enhancements	4
2.2	Backward compatibility of date formats in LIS communication	4
3	Bug fixes	5
3.1	Inability to select some items in correspondence tables	5
3.2	Correction of correspondence table errors for a deleted test	5
3.3	Improved handling of SIL connection errors	5

1 Object


This document provides a list of the modifications introduced in Version 16.02 of Valab® following Version 16.01.

2 Evolutions

2.1 LDAP connection enhancements

In the LDAP configuration window, it is now possible to specify multiple search paths in the dedicated field, by separating them with the '|' character.

Please note that the filter is shared across all search paths.

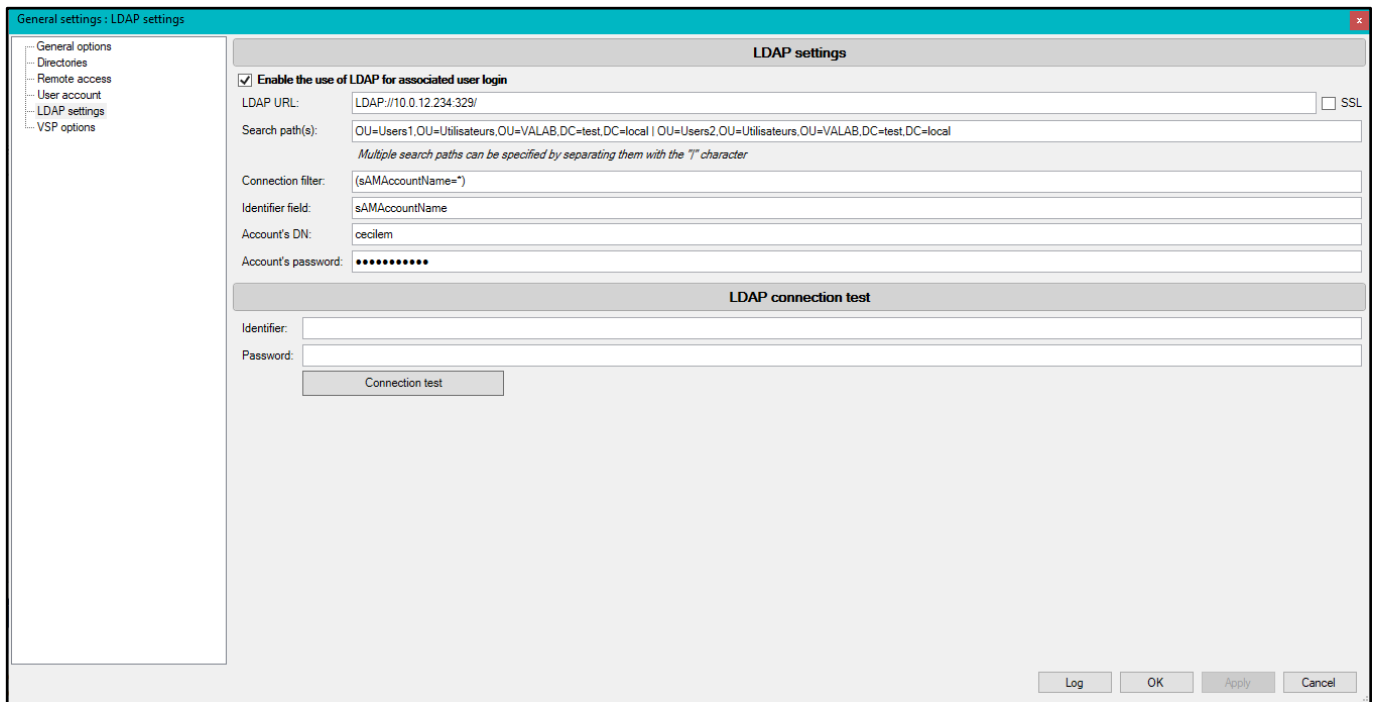

Specification: VLB-5155

2.2 Backward compatibility of date formats in LIS communication

An option has been added to maintain backward compatibility with some LIS systems using deprecated legacy date formats, with years represented by fewer than 4 digits or 6-digit dates (DDMMYY).

This option can be configured via an environment file and is disabled by default.

Specification: VLB-5163

3 Bug fixes

3.1 Inability to select some items in correspondence tables

The selection of items containing a dash or parentheses is now possible when managing correspondence tables.

Specification: VLB-5147

3.2 Correction of correspondence table errors for a deleted test

Error corrections for deleted symbolic test are no longer available in the correspondence table errors window.

Specification: VLB-5151

3.3 Improved handling of SIL connection errors

Low-level TCP/IP communication error handling with LIS has been improved to prevent potential blocking of the VALAB server.

Specification: VLB-5220